TERA ASOPS

High-Speed THz Time-Domain Spectrometer





With the ASOPS technique ultrafast data sampling is no longer limited by the natural restrictions of a mechanically moving optical delay unit. The novel TERA ASOPS high-speed THz time-domain spectrometer system is using this technique to temporally scan THz pulse traces at a revolutionary high rate. At the same time it is extending the detection window to nanoseconds pushing the spectral resolution into the region of hundreds of MHz.

TERA ASOPS is using two ultrafast lasers operating at a locked repetition rate with a tunable difference, providing the optical pulses for THz emission and detection. The lasers are featuring Menlo Systems' patented figure 9[®] mode locking technology for highest stability and reliability. The laser pulses are delivered via optical fiber to the THz antenna modules which are of the latest standard for high-power THz wave emission. A robust, plug-and-play off axis parabolic mirror terahertz optics ensures the highest bandwidth and long term stability of the THz signal for the TERA ASOPS. For seamless integration into existing experiments, the software features remote control of the spectrometer and high-speed data transfer over network. On request, additional laser output ports at 1560 nm and 780 nm can be configured.

PERFORMANCE DATA

THz pulse measured in low humidity atmosphere*

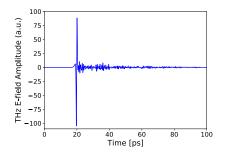

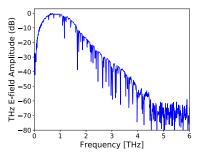

Calculated THz spectrum

*Settings: Averaging time: 500 s, Scanning rate: 100 Hz

MenioSystems

HIGHLIGHTS

- High-speed Data Acquisition
- Full Scanning Range
- Turnkey System Operation
- Intuitive, User-Friendly Software
- figure 9[®] Mode Locking
- Multi-Channel Option
- All-Fiber Flexibility
- Remote Control Engine

KEY SPECIFICATIONS

- >4.5 THz Bandwidth
- >70 dB Dynamic Range
- 10 ns Total Scan Range
- >600 Hz Scan Rate

APPLICATIONS

- Time Resolved THz Spectroscopy
- Chemical Fingerprinting
- THz Imaging
- THz Combs
- Stand-Off Detection

FEATURES

- Transmission and Reflection Geometry
- Fiber Coupled THz Path for Arrangement Outside the Spectrometer Housing
- Real-Time Measurements
- Remote Control via Network

OPTIONS

- Dual-Detection/Multi-Channel Suitable for parallel transmission and reflection measurements
- MULTIBRANCH Ask for Additional Optical Port Configuration
- TERA Image Automated XY Translation Stage for THz Imaging
- Reflection Head Compact Sensor Unit with Integrated THz optics
- THz Purge Box
 Enables Water Line Free THz
 Spectroscopy
- TeraLyzer Advanced Software for THz Data Analysis
- Polymer Lens Optics
- Custom Fiber Length

TERA ASOPS



High-Speed THz Time-Domain Spectrometer

SPECIFICATIONS

Spectral Range	>4.5 THz
Dynamic Range	>70 dB (typically 80 dB)
Total Scan Range	up to 10 ns (pulse-to-pulse distance)
Rapid Sampling Rate	>600 Hz
Laser Output Ports for THz	Two fiber coupled ports, 1560 nm, FC/APC, PM fiber, <90 fs after 2.5 m patch cord
System Repetition Rate	100 MHz (custom repetition rate available)

SYSTEM DIMENSIONS AND WEIGHT

Optomechanical Setup	900 x 600 x 200 mm ³ , 34 kg
THz Control Electronics	mounted in a 19" rack cabinet, 800 x 600 x 1800 mm ³ , 75 kg

SYSTEM COMPONENTS

Optical Breadboard	Two femtosecond laser sources	Two femtosecond laser sources C-Fiber*		
	Fiber coupled THz emitter and re	Fiber coupled THz emitter and receiver modules TERA15-FC*		
	THz TPX polymer lenses	THz TPX polymer lenses		
	THz off-axis parabolic mirrors op	THz off-axis parabolic mirrors optics		
TERA ASOPS Control Electronics	Laser control electronics	Laser control electronics		
	Synchronization electronics	Synchronization electronics		
	THz electronics	THz electronics		
	Data acquisition platform, 16 Bit	t, 105 MS/s		
	PC and software package for me	PC and software package for measurement and data analysis		
	22" screen, keyboard and mous	22" screen, keyboard and mouse		
	TCP socket remote control interf	face		
	.NET remote control interface	external analog/digital triggering		

*See product data sheet for detailed specifications.

REQUIREMENTS

Operating Voltage	110/115/230 VAC
Frequency	50 to 60 Hz
Cooling Requirements	no water cooling required
Operating Temperature	22 ± 5 °C

ORDERING INFORMATION	
Product Code	TERA ASOPS

Please call for pricing. Specifications are subject to change without notice. Custom modifications are available, please inquire.

MenioSystems

Menio Systems GmbH T+49 89 189 166 0 sales@meniosystems.com Menio Systems, Inc. T+1 973 300 4490 ussales@menlosystems.com **Thorlabs, Inc.** T+1 973 579 7227 sales@thorlabs.com



www.menlosystems.com

D-TERA_ASOPS-EN 21/06/21