

ORS

Ultrastable Laser System



MenloSystems

KEY SPECIFICATIONS

- Stability $<7 \times 10^{-16}$ at 1 s (ORS with ORS-FS-XTAL Option)
- Linewidth <1 Hz
- Output Power >10 mW
- Wavelength
650–1600 nm (ORS)
900–1600 nm (with ORS-FS-XTAL)

APPLICATIONS

- Optical Clocks
- Quantum Computing
- Ultra-low Noise Microwave Generation
- Frequency Comb Stabilization
- High Resolution Spectroscopy
- Laser Cooling and Trapping

OPTIONS

- **ORS-FS-XTAL:**
Crystalline Mirror Coating
- **ORS-EOM-SBL:**
EOM Sideband Lock
- **ORS-FNC:**
Fiber Noise Cancellation
- **ORS-MWL:**
Multiple-Wavelengths Mirror Coatings
- **ORS-SHG:**
Second Harmonic Generation
- **ORS-HP:**
High Output Power
- **ORS-DDRIF:**
Cavity De-drifting
- **ORS-DSRV:**
Digital Servo Interface for Stabilisation to External Atomic Transition

The Optical Reference System (ORS) is designed for uncompromised performance. It delivers ultra-narrow linewidth laser light with outstanding frequency stability.

The system's centerpiece is a high-finesse Fabry-Pérot cavity (cylindric spacer with a length of 12.1 cm) serving as a reference for a CW laser. The cavity is made out of ultra-low expansion glass (ULE) and is operated in vacuum at the point of zero thermal expansion. The reference cavity is actively decoupled from vibrations and acoustically isolated allowing for ultimate performance also in rough laboratory environments. An easy-to-use cavity-locking mechanism ensures portability without realignment of the optical paths.

The system is operated by the newest generation of Menlo Systems' proprietary SYNCRO controller—a modular electronics platform designed for versatility and intuitive use. It incorporates all required electronics, e.g., a low-noise laser driver and a very fast (analog) servo loop for laser frequency stabilization. The user controls all parameters using either the 7" front-panel touchscreen or a GUI on a remote PC. The onboard software ensures automatic cavity locking and system monitoring.

Ultrastable Laser System

SPECIFICATIONS


Wavelengths	650–1600 nm (ORS), 900–1600 nm (ORS with ORS-FS-XTAL)		
Stability (MADEV at 1 s, Linear Drift Removed)	<math> < 7 \times 10^{-16}</math> (ORS with FS-XTAL Option) <math> < 1 \times 10^{-15}</math> (ORS with ULE-IBS)		
Linewidth	<math> < 1</math> Hz		
Phase Noise (Laser Source Dependent)		ORS	ORS with ORS-FS-XTAL
	at 1 Hz	-8 dBc/Hz	-10 dBc/Hz
	at 10 Hz	-27 dBc/Hz	-30 dBc/Hz
	at 100 Hz	-47 dBc/Hz	-47 dBc/Hz
	at 1000 Hz	-70 dBc/Hz	-70 dBc/Hz
			For all models spurious signals <math> < -20</math> dBc
Output Power	>10 mW (FC/APC Connector), High Power on Request		
Cavity Spacer	12.1 cm, ULE, Cylindric Design		
Free Spectral Range	1.24 GHz		
Linear Drift Rate	approx. 150 mHz/s after 12 months		
Vibration Isolation Platform	Included		
Dimensions / Weight	590 x 800 x 1550 mm (29U) / 260 kg		

REQUIREMENTS


Operating Voltage	100 / 115 / 230 VAC
Line Frequency	50 to 60 Hz
Operating Temperature	22 ± 5 °C
Power Consumption	<math> < 150</math> W

STABILITY AND PHASE NOISE

ORS Fractional Frequency Stability
(ORS and ORS with ORS-FS-XTAL, 1542nm)

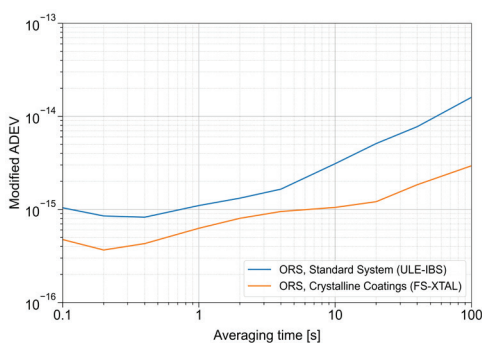

ORS with ORS-FS-XTAL Phase Noise (1542nm)

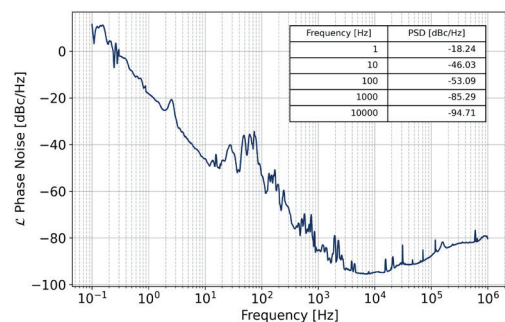

ORDERING INFORMATION

Product Code	ORS
--------------	-----

Please call for pricing. Specifications are subject to change without notice. Custom modifications are available, please inquire.



Invisible laser radiation
avoid exposure to beam
Class 3b laser

Menlo Systems GmbH
T+49 89 189 166 0
sales@menlosystems.com

Menlo Systems US
T+1 303 635 6406
ussales@menlosystems.com

Menlo Systems Japan
T+81 907 409 20 21
jpsales@menlosystems.com

Menlo Systems China
T+86 21 6071 1678
chinasales@menlosystems.com

